

PROTOCOL K-300i - DILUTION DIRECTIONS

ATTENTION

K-300i IS A CONCENTRATED PRODUCT AND MUST BE DILUTED WITH WATER. PLEASE FOLLOW THE DIRECTIONS BELOW.

EACH BOX CONTAINS

- ✓ Four 4L jugs of CONCENTRATED K-300i
- ✓ One 4L calibrated pump
- ✓ Envelope containing (ready to use) labels for 1L bottles of diluted K-300i
- ✓ Dilution directions
- ✓ Product Information Sheet

DIRECTIONS

1. Assemble the pump by inserting flat end of the straw into the bottom of the pump.
2. Open a 4L jug of CONCENTRATED K-300i and attach pump by inserting and screwing down tightly to the top of the jug.
3. Unlock pump by pushing down and twisting pump to the left.
4. Add 1-3 pumps of CONCENTRATED K-300i to an empty 1L bottle (not provided) and fill the rest with water.
5. Attach provided label to the 1L bottle and attach a spray nozzle (not provided).
6. You are ready to clean! See product information sheet for application instructions.